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Diamond Products

The evolution of diamond products:

Diamond Paste was the first diamond product to be used instead of alumina for polishing of metallographic samples. A small amount of diamond paste is distributed manually on the polishing cloth before preparation. During the preparation process only lubricant is added to achieve the correct lubrication and cooling action. Diamond paste cannot be added during the preparation process, it cannot be dosed automatically and it cannot be used on rigid grinding discs.

When Diamond Suspension was developed, precise automatic dosing was made possible by using peristaltic pumps and reproducible preparation methods with well-defined dosing levels were a reality. However, as from the beginning, diamond and lubricant were still applied individually.

The next logical step was to combine suspension and lubricant into one liquid, the 2-in-1 Diamond Suspension. These 2-in-1 suspensions guarantee the correct suspension/lubricant ratio at all times and combine the best possible preparation results with utmost ease of use.

In between the development of the different suspensions the Diamond Stick was introduced as a substitute of the diamond paste. The diamond stick is easier to use than diamond paste, and is, because of the very high diamond concentration, recommended for priming new, unused polishing cloth. It can, contrary to Diamond Paste, also be added during preparation.

To cater for all the different requests and to satisfy our customers' wishes, we have developed the widest selection of high quality diamond products.

Most of our diamond products are available in two different versions; with either polycrystalline or monocrystalline diamonds.

The first diamonds to be used for metallographic polishing were natural, monocrystalline diamonds, as used in the production of jewellery. Since then the demand for diamonds for polishing applications has increased dramatically and diamonds are now produced artificially.

Artificial monocrystalline diamonds have replaced natural diamonds in most metallographic applications because they have the same properties. For the highest requirements and the best preparation results, polycrystalline diamonds are used.

Polycrystalline diamonds break down over time, continuously producing small abrasive edges that provide high material removal while producing only shallow scratches. Monocrystalline diamonds will get more and more round over time, losing their cutting ability.

2-in-1 Suspensions	DiaUltra	DiaMaxx		DiaDoublo	
	Mono	Poly	Mono	Poly	Mono
Ease of use	✓✓✓✓✓	✓✓✓✓✓	✓✓✓✓✓	✓✓✓✓✓	✓✓✓✓✓
Diamond concentration	+++	+++	+++	++	++
Removal rate	↑↑↑↑↑	↑↑↑↑	↑↑↑	↑↑↑	↑↑
For water-free preparation	÷	÷	÷	÷	÷

Suspensions	Aka-Poly			Aka-Mono	
	Aka-Poly+	Aka-Poly	Aka-Poly WF	Aka-Mono+	Aka-Mono
Ease of use	✓✓✓✓	✓✓✓✓	✓✓✓✓	✓✓✓✓	✓✓✓✓
Diamond concentration	+++	++	++	+++	++
Removal rate	↑↑↑↑↑	↑↑↑	↑↑↑	↑↑↑↑	↑↑↑
For water-free preparation	÷	÷	✓	÷	÷

Other	Diamond Stick		Diamond Paste	
	Poly	Mono	Poly	Mono
Ease of use	✓✓	✓✓	✓	✓
Diamond concentration	++++	++++	+++++	+++++
Removal rate	↑↑↑	↑↑↑	↑↑↑	↑↑↑
For water-free preparation	✓	✓	✓	✓